


Twin-Strep-tag Capture Kit

with red dye **Ra**

Dynamic Biosensors GmbH & Inc.

HK-SXT-1 v6.1

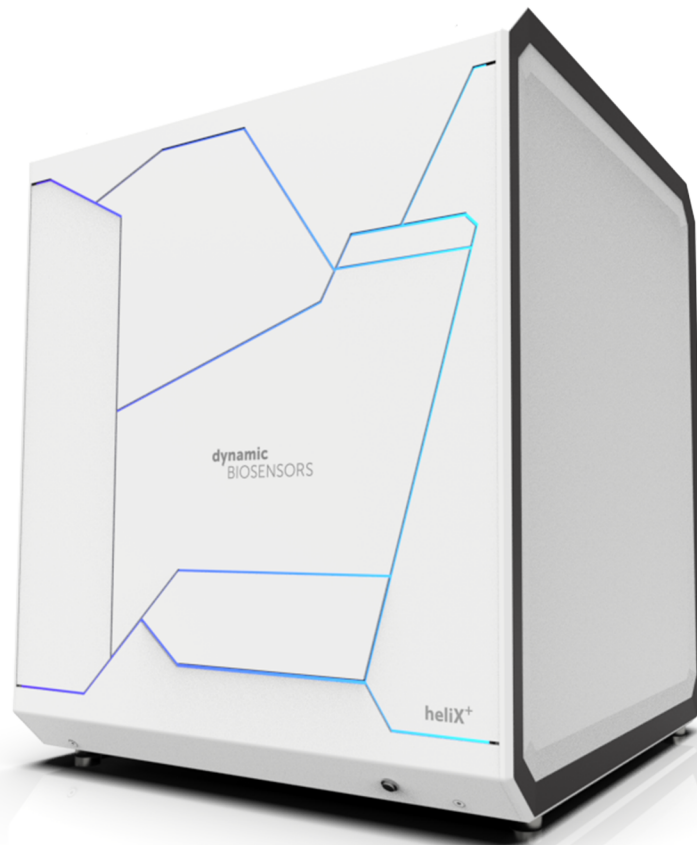


Key Features

- This kit is designed for capture of **Twin-Strep-tagged molecules** using **Strep-Tactin[®]XT (SXT)**.
- Compatible with **heliX[®] Adapter Chip**.
- Includes **Adapter strands** and **Ligand strand** modified with Strep-Tactin[®]XT for **20 regenerations**.
- For functionalization of **Spot 1** and **Spot 2**.
- **Adapter strands 1** and **2** carry a moderately hydrophilic red dye (**Ra**) with a single positive net charge.

heliX[®] Adapter Chip Overview

2 spots with 2 different anchor sequences for DNA-encoded addressing. Spot 1 is functionalized with the capture molecule while Spot 2 is used as real-time reference.

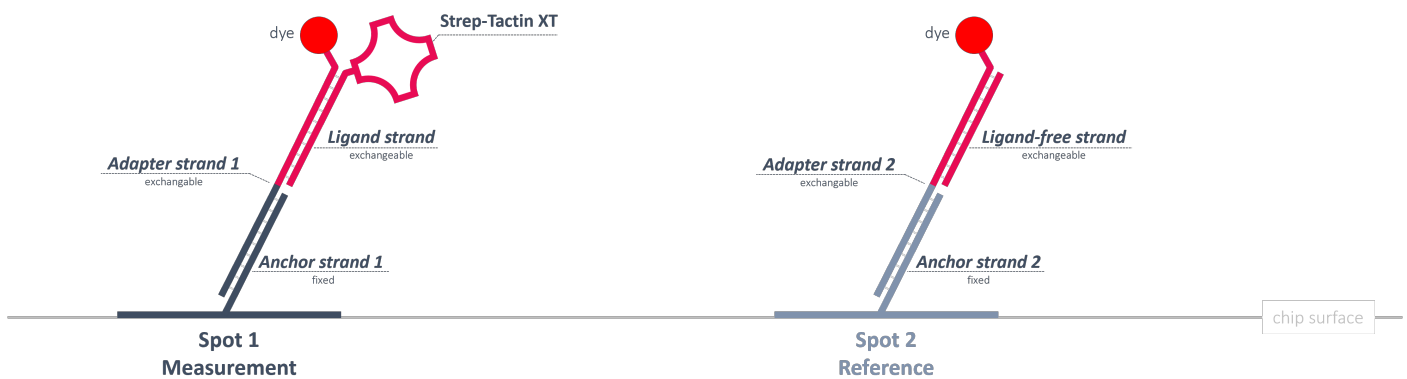

Product Description

Order Number: **HK-SXT-1**


Table 1. Contents and Storage Information

Material	Cap	Concentration	Amount	Buffer	Storage
SXT^[1] - Ligand strand	Green	500 nM	2 x 100 µL	TE40 ^[2]	2-8°C
Adapter strand 1 - Ra	Black	400 nM	2 x 100 µL	TE40 ^[2]	-20°C
Adapter strand 2 - Ra - lfs	White	200/250 nM	2 x 200 µL	TE40 ^[2]	-20°C

For research use only.

This product has a limited shelf life, please see expiry date on label. After preparation of ready to use solution the expiry date is **3 months**.

Workflow of a heliX[®] SXT - capture assay

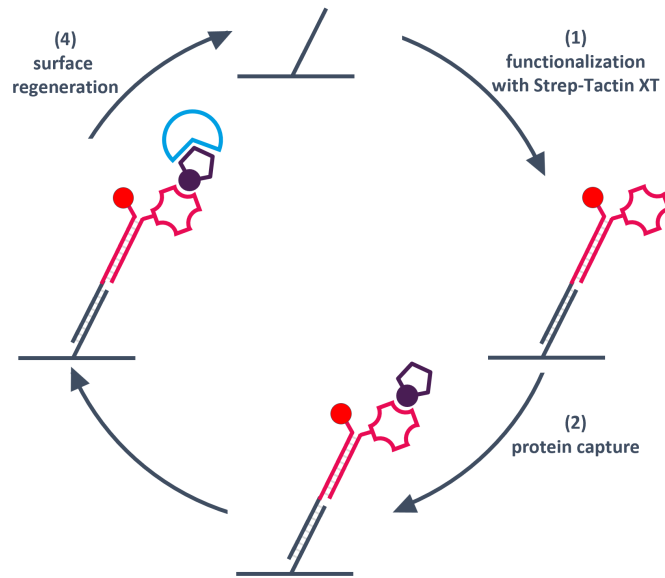

1. The anchor strand (ssDNA) immobilized to the surface of the **heliX[®] Adapter Chip** is hybridized with complementary DNA strands modified with **Strep-Tactin[®] XT**.
2. The **twin-strep-tagged ligand** of interest is captured on the surface during the measurement run.
3. Measurement of the analyte binding kinetics or conformational change upon analyte binding.
4. Surface regeneration by injection of a high pH solution. Chip surface goes back to the original state. This step can be followed by a new hybridization of fresh ligand with Strep-Tactin[®] XT.

Preparation

1. Mix 100 μ L **SXT - Ligand strand** with 100 μ L **Adapter strand 1 - Ra**.
2. Incubate the solution of step 1 at **RT** at **600 rpm** for **30 min** to ensure complete hybridization.
3. Mix 200 μ L **Adapter strand 2 - Ra - lfs** to the sample after step 2.

The solution (400 μ L in total) is ready to use for a biochip functionalization.

Please aliquot and store the ready to use solution at 2-8°C. **Use up within 3 months.**

The kit contains material for the preparation of two separate ready to use solutions with 400 μ L each.

Assay Setup in heliOS

Go to **heliOS** > create a **New Assay Workflow** > add **Custom Assay** > load **Capture with Kinetics** > modify the parameters based on your needs and run the assay.

Suggested assay parameters (e.g., flow rate, time, LED power, etc.) are within the **heliOS** assay.

For further questions, please contact the support team at support@dynamic-biosensors.com

Useful Order Numbers

Table 2. Order Numbers

Product Name	Comment	Order Number
heliX[®] Adapter Chip	Chip with 2 detection spots	ADP-48-2-0
10x Passivation solution	For passivation of chip surface	SOL-PAS-1-5
Regeneration solution	For regeneration of chip surface	SOL-REG-1-5

Contact

Dynamic Biosensors GmbH

Perchtinger Str. 8/10
81379 Munich
Germany

Dynamic Biosensors, Inc.

300 Trade Center, Suite 1400
Woburn, MA 01801
USA

Order Information order@dynamic-biosensors.com

Technical Support support@dynamic-biosensors.com

www.dynamic-biosensors.com

Instruments and chips are engineered and manufactured in Germany.

©2024 Dynamic Biosensors GmbH | Dynamic Biosensors, Inc. All rights reserved.

[1] Strep-Tactin®XT is a product of IBA GmbH

[2] TE40: 10 mM Tris, 40 mM NaCl, 0.05 % Tween20, 50 µM EDTA, 50 µM EGTA