


switchSENSE[®] Biochip MPC2-96-2-G1R1-S

Product Description

Product Code	MPC2-96-2-G1R1-S
Suitable for	switchSENSE [®] DRX ^{red} & DRX ²
Storage	Store at 2-8 °C, dry in the dark
Layout	6 double reference spots (nanolever A/B)
Grade	Standard

Bi-functional chips offer a very flexible and powerful platform to study bi-functional antibodies, e. g. to discriminate between complex binding modes (monovalent versus bivalent or bispecific).

Biochip Design

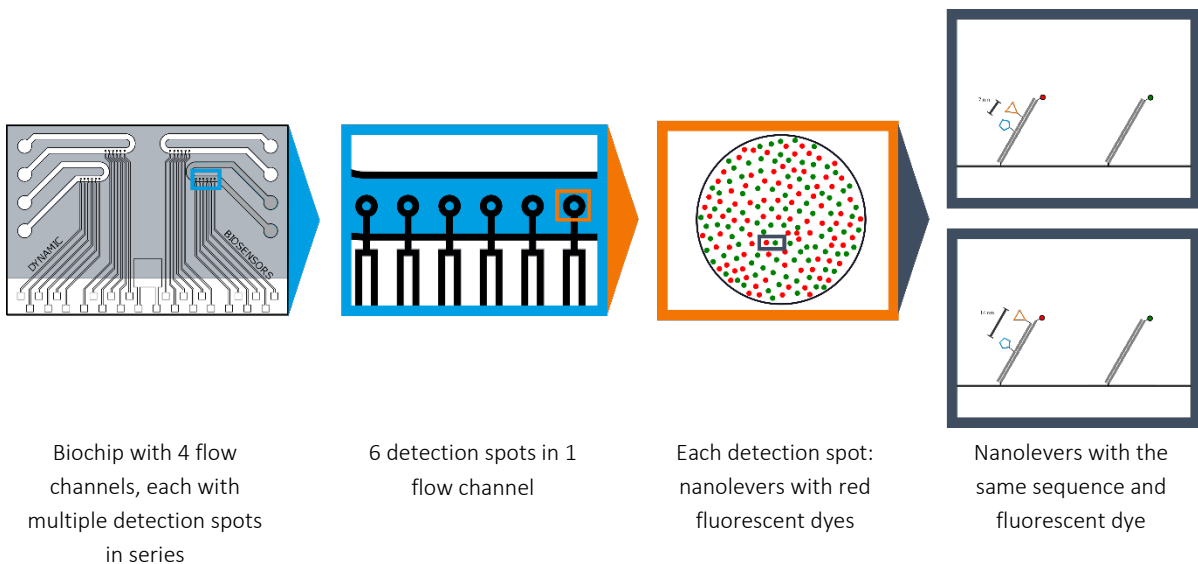


The detection spots are functionalized with two different nanolevers, NL-A96 and NL-B96. The 5'-end of the DNA is fixed on the surface while the 3'-end carries a fluorescent dye, either a green one (G1) on NL-A96 or a red one (R1) on NL-B96.

For functionalization either CK-NH2-5-B96 or CK-NH2-6-B96 are available. The first kit provides the possibility to bind two antigens in a distance of 7 nm on the nanolever B96, the later features a distance of 14 nm for the two antigens.

Flow channel with 6 double (A, B) spots (real-time on-spot referencing or sophisticated assays)

BIF2 biochips carry a mixture of two nanolever populations (red & green) on the same sensor spot. Use one nanolever as the target and the other as the control for absolute data confidence. Beyond that one can use this set-up to investigate more complex and challenging issues.

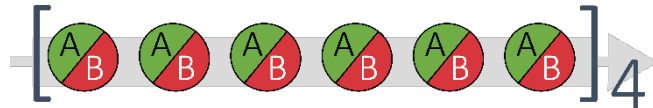

For more information about the DNA sequences, please contact us at info@dynamic-biosensors.com.

Application Areas

<i>Binding Kinetics*</i>	k_{on}, k_{off}, K_D
<i>Binding Affinity*</i>	K_D , titration curve, n (Hill coefficient)
<i>Protein Diameter*</i>	D_H (Hydrodynamic diameter)
<i>Conformational Change*</i>	ΔD_H (relative change of hydrodynamic diameter)
<i>Melting & Thermodynamics*</i>	$T_M, \Delta G, \Delta H, \Delta S$
<i>Multimers & Aggregation*</i>	Monomer-dimer discrimination, aggregation
<i>Nuclease & Polymerase Activity*</i>	$k_{cat}, K_M, T_{activate}$
Bispecific Binders & Avidity	Binding affinity/avidity/kinetics

* for this application, other biochips (MPC or ENZ) are recommended

Biochip Handling

The biochip is ready to use. For research only.

Avoid touching or picking up the biochip with your hands as this may cause bilayer electrostatic discharge. Instead, always use the vacuum tweezers provided with the instrument when handling the biochips.

After installing the biochip in the DRX instrument and before starting an experiment be sure that the selected channel is passivated with passivation solution (Order No. SOL-PAS-1-5).

More information about biochip storage and handling can be found in the document “switchSENSE Biochips Storage Information”.

Grade I Standard

Standard grade chips are batch tested which means one out of a batch undergoes a comprehensive quality control – proper switching dynamics, calibration (voltage-response of the DNA layer) and fluorescence levels. In addition each biochip is tested for proper fluorescence levels on each electrode. Due to the production process, not all detection spots fulfill our high quality standards. Each standard grade biochip is certified as having 20 or more active detection spots.

Compatible Functionalization Kits

All compatible functionalization kits for this biochip are listed on our website:

www.dynamic-biosensors.com/reagents/

Contact

Dynamic Biosensors GmbH

Lochhamerstr. 15
82152 Martinsried/Planegg
Germany

Order Information

Phone: +49 89 89 74 544 0
order@dynamic-biosensors.com

Dynamic Biosensors Inc.

9705 Carroll Centre Road, Suite 100
San Diego, CA 92126
USA

Technical Support

Phone: +49 89 89 74 544 66
support@dynamic-biosensors.com